DH Kids DMS1

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Spis treści

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
ENGLISH

Table of contents

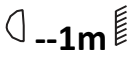Safety rules.....	9
Maintenance	10
Operation	10
Information about used electrical and electronic equipment	14

Safety rules

	Please read the user manual in detail before use. Please also keep it for future reference. The manual contains rules for the safe use of the device.
	After receiving the product, please unpack and check whether it is complete and whether there is any damage caused by transportation. If there is damage caused by transportation, please do not use this product and contact the dealer or manufacturer as soon as possible.
	WARNING! THE DEVICE MUST NOT BE DISPOSED OF WITH HOUSEHOLD WASTE. This symbol indicates that this product must not be disposed of with household waste, according to the EU and your national law. In order to prevent potential damage to the environment or health, the used product must be recycled. In accordance with current legislation, unusable electrical and electronic devices must be collected separately at the designated facilities for recycling, acting on the basis of applicable environmental standards.
	The product described in this manual comply with European directives and it is therefore CE marked.
	Keep the device away from children and unqualified persons. The manufacturer is not liable for damage caused by improper use.
	Before use, make sure that the housing is not damaged.
	Ensure a minimum distance of 1 m between the appliance and flammable materials.
	The device operates on the power supply marked on the housing - do not connect to a power supply with different parameters. Disconnect the product from the power supply by pulling on the power plug - do not pull directly on the power cord! Disconnect the device from the power supply before removing the housing or performing maintenance. If you find the power cord is damaged - do not use it!
	Indoor device, use in a well-ventilated room, keep away from moisture. The distance between the device and a wall should be kept at least 50 cm. Do not expose the product to direct sunlight or lighting devices. Do not store or use the device near a fire source!
	Do not look directly into the light source - the light may cause an epileptic attack in light-sensitive people or people with epilepsy. Compliance with this rule is mandatory.
	Do not install the device on a surface subjected to vibration. Do not expose the device to high temperatures!
	In the event of a malfunction, disconnect the device from the power supply immediately!


Maintenance

1. The device may only be used by qualified personnel, damage caused by improper use or attempted repair is not covered by the warranty. There are no service parts inside the packaging, repairs may only be carried out by authorized service center.
2. Do not allow light source to come into contact with oil, grease or any similar liquid. The light source should not come into direct contact with the skin!
3. Regular cleaning allows for long-term use and maintains proper quality of the device operation. Use a soft cloth to clean the device.

Operation

1. Connect the device to the power source, observing the safety rules of use, and fully charge it. Start the fully charged device with the power button.
2. The device shall be operated manually only.
3. When you have finished using the device, turn it off with the power button and disconnect it from the power source, observing the safety rules for use.

Structure

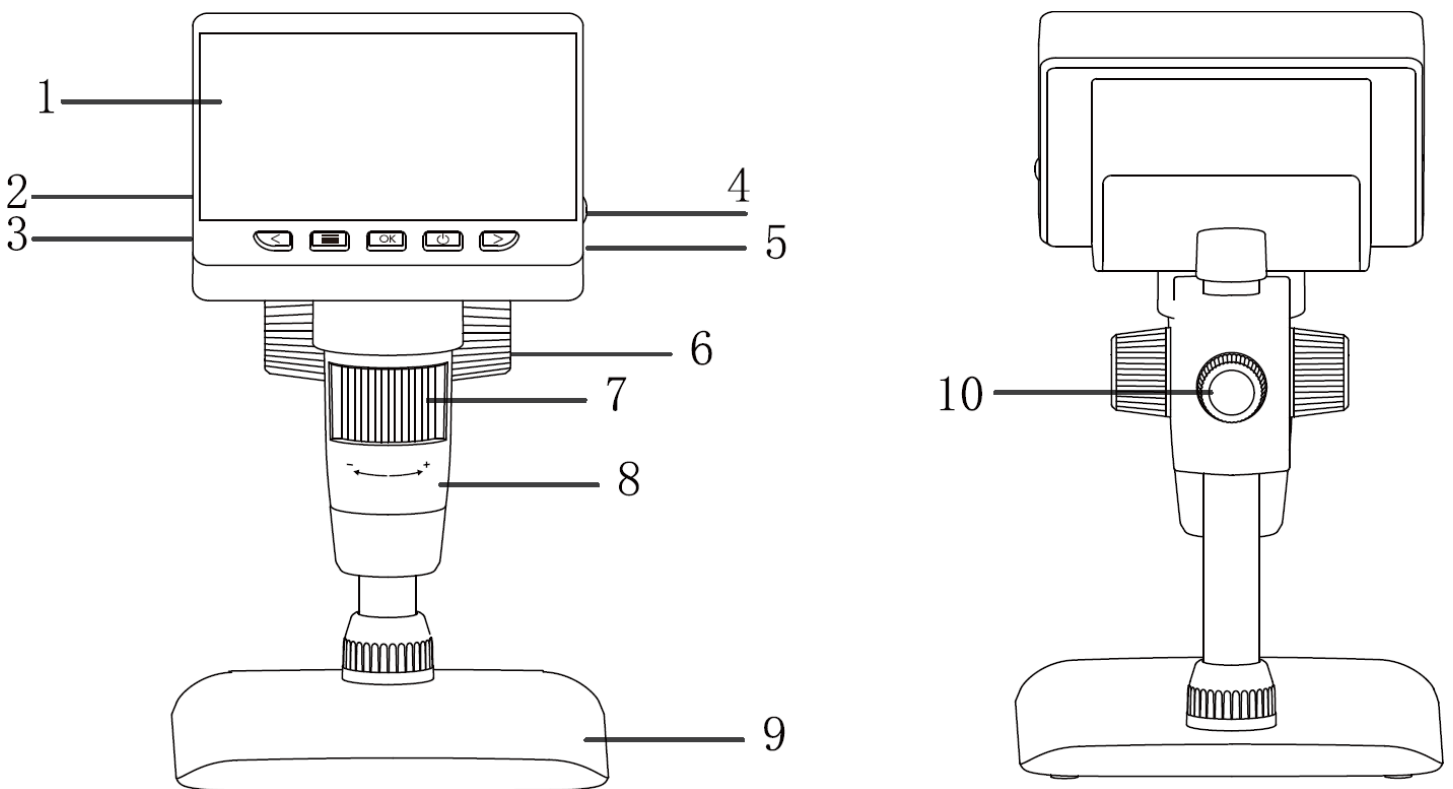1. 4.3-inch display.
2. MicroSD card port.
3. USB power input.
4. LED light source adjustment knob.
5. Reset hole.

6. Knob A.
7. Focus adjustment knob (knob C).
8. Tube.
9. Base.
10. Knob B.

Control panel and basic functions

1. Zoom out/move up/left button - in photo mode, a short press zooms out and a long press rotates; in other modes, a press moves up or left.
2. Menu/return button - short press to go to the advanced settings screen or return to the photo preview screen; in the advanced settings screen, you can view or set the interface language / date and time / screen saver / light source frequency / format / default settings / version.
3. OK button - in Photo mode, short press to take a photo; long press to enter video recording mode - short press again to start recording and short press to stop recording.
4. Power button - short press to enter and exit file playback mode; long press to turn the device on or off.
5. Zoom in/move down/right button - in photo mode, short press to zoom in; in other modes, press to move down or right.

MicroSD card port - supports card size up to 32 GB with FAT32 format; microSD card is not included.

USB power input - DC 5V 1A input to provide power or data transfer.


Focus adjustment knob - turn the knob to control the focus of the image.

LED light source adjustment knob - turn the knob to control the brightness of the light source.

Reset hole - pressing the button at the bottom of the hole allows you to reset the device.

Use

1. Insert the microSD card (optional) into the card port. Gently push in until the card is locked. Note: the microSD card must be inserted to record and store photos and videos.
2. Press and hold the power button to turn on the device.
3. Place the object to be observed on the stage, turn knobs A and B to bring the object closer to the microscope tube.
4. Turn knob C to the left or right to adjust the focus precisely.
5. When the desired focus is in place, turn knob C to lock the stage.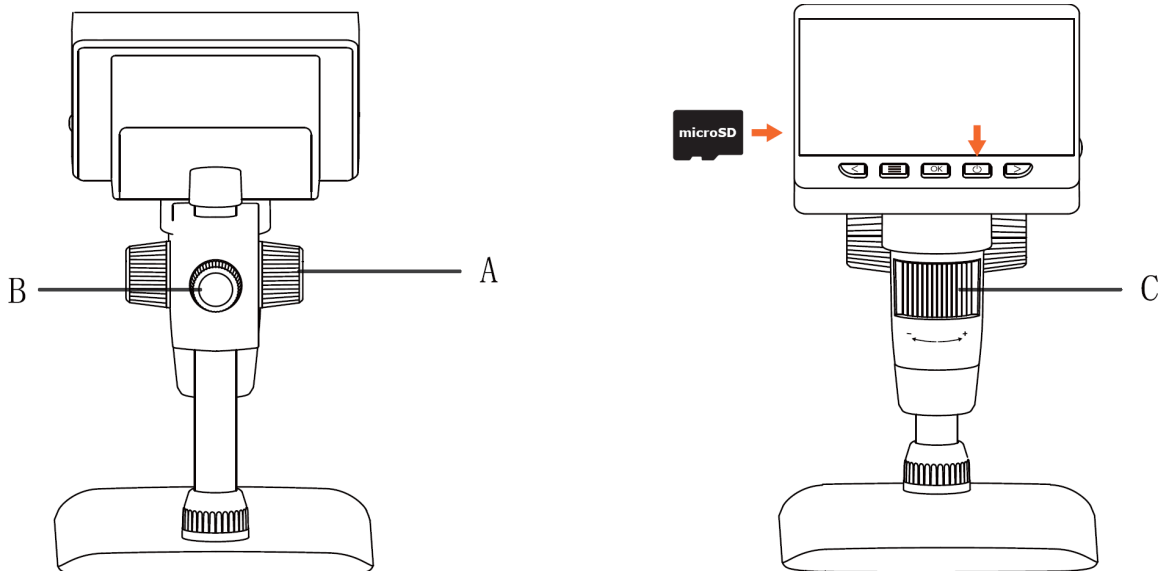

Operation modes



The device is equipped with two operation modes:

- memory mode: connect to the computer to load, save and format the microSD card; connect the microscope to the computer using the USB data cable, select the storage mode, and then click the OK button to display the saved photos and videos on the computer (the device control panel buttons will not be available in storage mode - only the power button and the LED light source adjustment knob will remain active);
- PC camera mode: connect to a computer and use the device with computer software; in PC camera mode, the device menu button will not be active, the device will display a blue screen in PC camera mode (which is a normal phenomenon, not a problem with the device), and if there is no image in the software, repeat connecting and disconnecting the USB data cable.

PC camera mode


Windows

1. For Windows 10/11 systems, it is recommended to use the Camera app, but you should disable the default system camera and grant appropriate camera privacy permissions. For Windows 7/8 systems, it is recommended to download and install AMCAP software.
2. Plug the USB cable into the USB port on the back of the computer case. Press and hold the power button to turn on the device and select PC camera mode.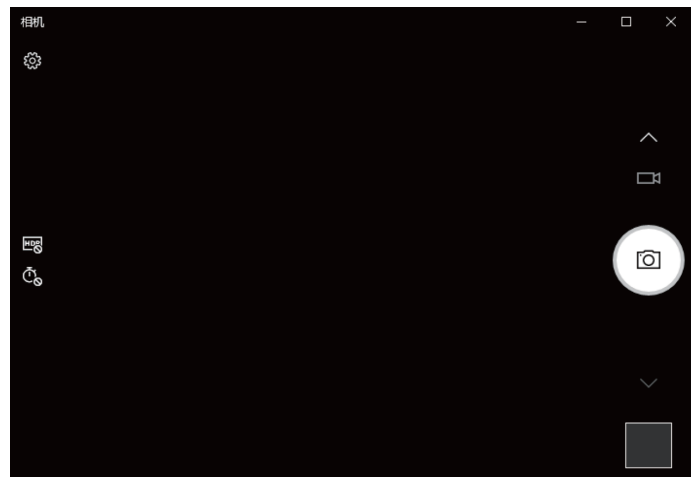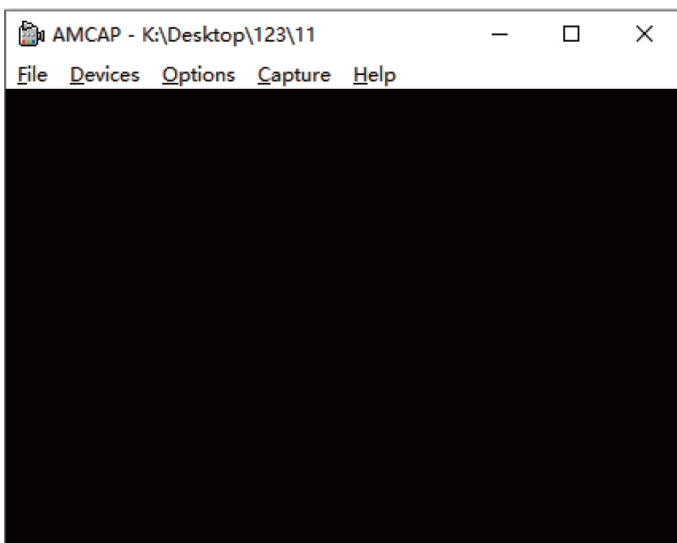


MacOS

1. For MacOS 11 and above, it is recommended to use Photo Booth software.
2. Plug the USB cable into the USB port on the back of the computer case. Press and hold the power button to turn on the device and select PC camera mode. Open the Photo Booth software.
3. Select "Camera" and then "USB CAMERA."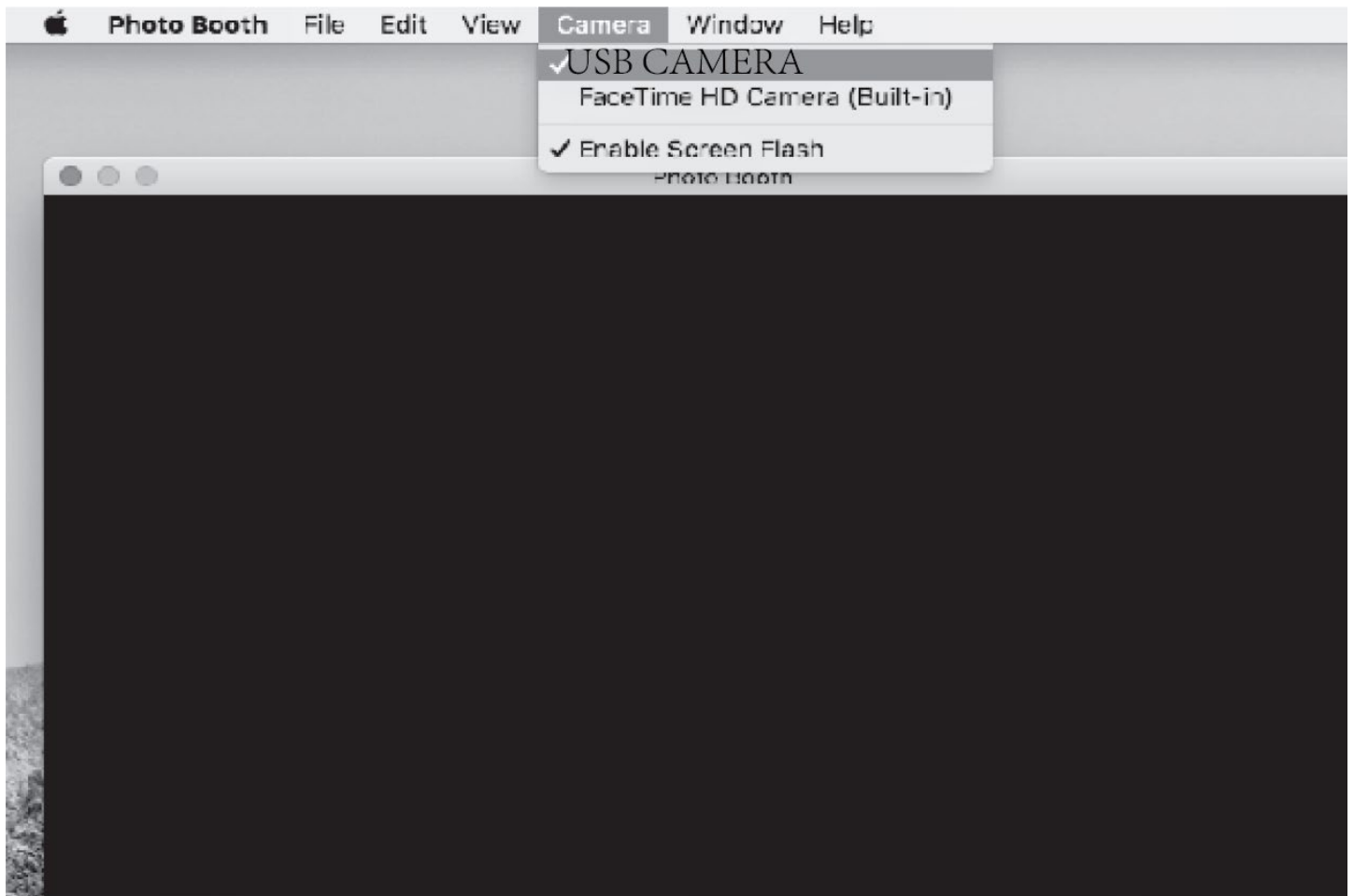

Charging

1. Use the 5V 1A charger adapter to charge the device.
2. During charging, an icon will appear on the microscope display. When charging is complete, the battery charge indicator looks as shown below.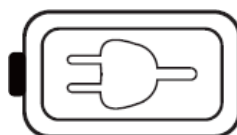

3. When the device is fully charged, the microscope can be used continuously for up to about 6 hours.
4. After the device is fully discharged, charge it with a 5V/1A charger for at least 1 hour before you can use it normally.
5. Do not charge the device with a computer. The device will automatically go into PC camera mode.

Information about used electrical and electronic equipment

The main goal of European and national law regulations is to reduce the amount of waste produced from used electrical and electronic equipment, to ensure an appropriate level of collection, recovery and recycling of used equipment, and to increase public awareness of its harmfulness to the environment, at each stage of use of electrical and electronic equipment. Therefore, it should be pointed out that households play a key role in contributing to reuse and recovery, including recycling of used equipment. The user of electrical and electronic equipment - intended for households - is obliged to return it to authorized collector after its use. However, it should be remembered that products classified as electrical or electronic equipment should be disposed of at authorized collection points.